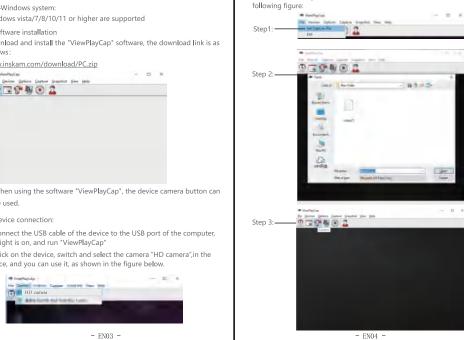
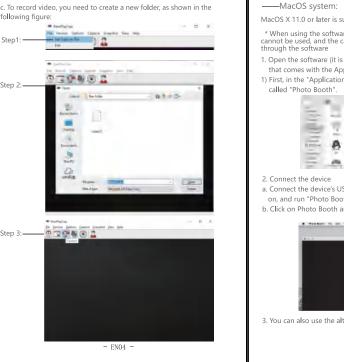
Product description ——Computer users Preface ----Windows system: Thank you for choosing this equipment. The usage of this product is Windows vista/7/8/10/11 or higher are supported complicated, please read this manual carefully before using to understand 1, Software installation CONTENTS the operation of microscope. Download and install the "ViewPlayCap" software, the download link is as follows: Important Notes 1. Camera button www.inskam.com/download/PC.zip 1. Please open the dust cover in front of the lens when using it for the first - D X English..... —2 2. Light adjustment key Die Gerices Options Capture Graphot 2. The camera button on the product can only be used on the software 3. Focusing tube "ViewPlayCap", other software cannot be used. Hello —3 4. Spotlight Cup 3. This equipment is a combination of digital microscope and optical Chinese..... ..08-14 microscope. The specific magnification effect is subject to the actual 5. Dust cover shooting picture. 4. The best focal length of the microscope is 0-40mm. The image is not clear beyond 40mm, you need to adjust the focus by adjusting the focus wheel to see the object clearly. Product introduction * When using the software "ViewPlayCap", the device camera button can Thank you for purchasing our WiFi electron microscope, which can be easily be used. used in different areas including: 2. Device connection: Textile industry, used for textile inspection a. Connect the USB cable of the device to the USB port of the computer, User Manual the light is on, and run "ViewPlavCap" 3. PCB, precision machinery testing b. Click on the device, switch and select the camera "HD camera",in the 4. Educational purpose device, and you can use it, as shown in the figure below. 5. Hair check Thank you for choosing 6. Skin examination digital microscope 7. Microbial observation 8. Inspection of jewelry and coins (collection) 9. Visual AIDS

Frequently Asked Questions MacOS X 11.0 or later is supported. * When using the software "Photo Booth", the camera button of the device cannot be used, and the camera and video function can only be realized through the software A1: Support Windows Vista/7/8/10/11, Mac OS X 11.0+, Linux/Chrome OS is not supported. 1. Open the software (it is recommended to use the "Photo Booth" software

that comes with the Apple computer). 1) First, in the "Applications" directory of the Finder window, find a software

called "Photo Booth". BATT - - - -2. Connect the device

a. Connect the device's USB cable to the computer's USB port, the light is on, and run "Photo Booth". b. Click on Photo Booth and select the camera "HD camera",

President Sk. Str. View | _____ Stroke Skip
HD camera

3. You can also use the alternative software "Quick Time Player".

- EN05 -

- Q1: What systems is this device compatible with?
- O2: Liust got it and connected it to my computer, it looks like the camera
- button doesn't work? A2: The camera button on the product can only be used on the software
- Q3: Why is the image very blurry when viewing?

"ViewPlayCap", other software cannot be used.

- A3: 1. Have you adjusted the microscope to the best observation distance You can adjust the focus by slowly rotating the focusing wheel left and right to achieve the clearest state.
- 2. Whether to open the dust cover in front of the lens.
- Q4: What is the reason for the unmoving black patches displayed on the screen when using the device to watch?
- A4: Due to some reasons, dust has entered. You can try to tap the produc to see if the dust can be knocked out. If not, return it to the factory for
- and opening the software, there is no image. A5: 1. Reconnect the microscope and PC side.

Q5: The microscope is connected to the PC, after successfully downloading

2. Connect to the USB port on the back of the computer host.

- EN06 -

Q6: After the device is successfully connected, what is the reason why the LED light does not light up and the display has no image?

- A6: 1. Confirm whether the light is turned off.
- Reconnect the microscope and computer.
- O7: After the device is successfully connected, what is the reason why the LED light is on and the display has no image?
- A7: The software used is not compatible with this computer system. Alternative software is optional.
- Q8: Can the device record sound?
- A8: No, it cannot record sound.

Technical Parameters:

Pixel	0.3 million pixels
Magnification	50X-500X/50X-1000X/50X-1600X
Picture resolution	640*480P
Focal distance	Manual focus(0~40mm)
Picture format	JPG
Video format	AVI
Dynamic frame rate	20-30f/s
USB interface type	USB 3.0/2.0/1.0
Supportive system	Windows vista/7/8/10/11+ , Mac OS X 11.0+

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